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QMOUNT

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UV 埋込機 QMOUNT  
**SUPERIORITY IN DETAIL**

### 1. Robust housing

- | Sturdy powder-coated aluminum construction
- | Easy to clean
- | Integrated ventilation

### 2. Large scratch-resistant glass plate

- | Allows up to 12 samples simultaneously
- | 外形寸法: 200 mm x 260 mm]

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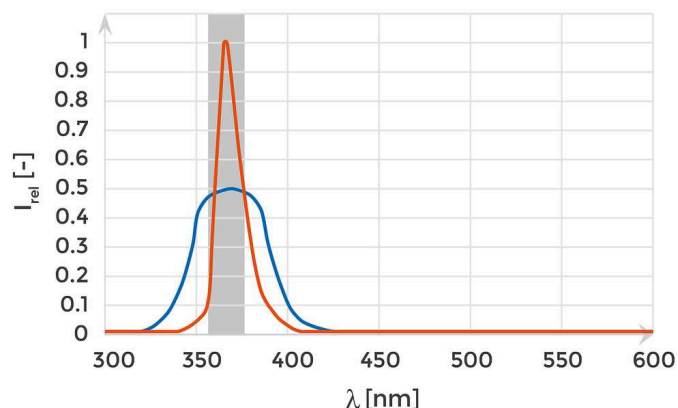
To minimize loss of electric power the Qmount is equipped with specialized UV LEDs. These show a very slim emission spectrum, which maximum superimposes with the absorption maximum of the KEM 50 UV initiator. This enables shorter curing times with a low electric consumption.

UV initiators get the necessary activation energy for the initiation reaction by absorbing UV radiation. Every molecule has a specific absorption spectrum, which shows certain wavelength areas in which the absorption coefficient of the molecule is particularly large. Outside of the so called absorption maxima a large portion of the emitted radiation is lost. Standard UV LEDs usually emit a significant amount of radiation besides the wavelength intervals in which the initiator has a good absorption. This radiation is converted to heat.

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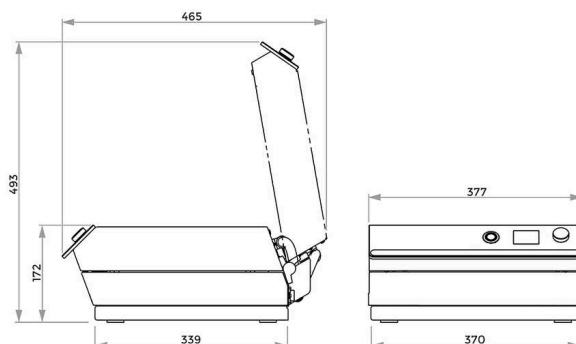
[www.qatm.com/qmount](http://www.qatm.com/qmount)

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